## Botanical Reference Material Product Grades

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### Reagent Grade Botanical Reference Materials (RBRMs™)

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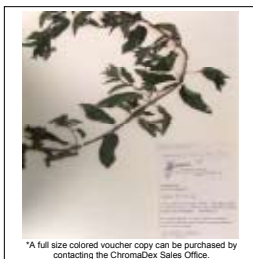
# Example: VBRM™ Certificate of Analysis

ChromaDex certificate of analysis

PRODUCT	<i>Mentha piperita</i> (Peppermint) Leaf
PART NUMBER	00030777
REFERENCE TYPE	Biomass Reference Material (V)*
LOT NUMBER	
CDXA NUMBER	
LATIN NAME	<i>Mentha piperita</i> L. [Lamiaceae]
COMMON NAME	Peppermint
PLANT PART	Leaf
VOUCHER NUMBER	
REPORT NUMBER	
DATE OF SAMPLE	14 July 2003
DATE OF REPORT	22 May 2007

\*A full size colored voucher copy can be purchased by contacting the ChromaDex Sales Office.

ChromaDex certificate of analysis

## MACROSCOPY

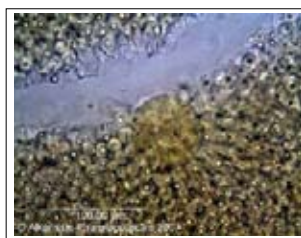
## ANALYTICAL CONDITIONS

TEST	METHOD	RESULT
Macroscopy	Digital Image	Green and beige fragments of leaf
Microscopy	Digital Image (400X)	(1) Surface of the leaf shows glandular trichomes with 8 secretory cells (2) Fragment of large covering trichomes with striated cuticle
HP-TLC	CDXA-TLCM-014-00	Conforms (See Below)
Appearance	Visual (Milled)	Green fluffy powder

## STORAGE CONDITIONS

STORAGE	20-30 °C; Dry storage area; Insect free; Volatile free
EXPIRATION DATE	05/2012 under the above conditions.

## MICROSCOPY


ChromaDex certificate of analysis

## HP-TLC CONDITIONS

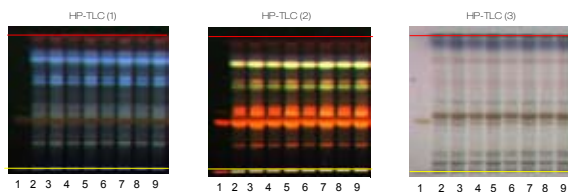STATIONARY PHASE	Silica gel 60, F <sub>254</sub> , 10 x 10 cm HP-TLC plates	
SAMPLE PREPARATION	0.5 g + 5 ml CH <sub>3</sub> OH sonicated in hot water ~ 10 minutes, heat in oven @ 65° C ~ 30 min.	
MOBILE PHASE	Ethyl acetate/ Acetic acid/ Formic acid/ Water	[10/1.1/1.1/2.6]
CHAMBER TEMPERATURE	Ambient	
DETECTION	(1) UV 365 nm (2) Natural Products spray reagent → UV 365 nm (3) Vanillin/H <sub>2</sub> SO <sub>4</sub> spray reagent → visible light	

## HP-TLC CHROMATOGRAMS

YELLOW LINE = SAMPLE ORIGIN  
RED LINE = SOLVENT FRONT @ 60MM

## HP-TLC APPLICATIONS

Lane	ID	Lane	ID
1	Rutin	6	<i>Mentha piperita</i> Leaf VBRM
2	<i>Mentha piperita</i> Leaf VBRM	7	<i>Mentha piperita</i> Leaf VBRM
3	<i>Mentha piperita</i> Leaf VBRM	8	<i>Mentha piperita</i> Leaf VBRM
4	<i>Mentha piperita</i> Leaf VBRM	9	<i>Mentha piperita</i> Leaf VBRM
5	<i>Mentha piperita</i> Leaf VBRM		

## General Overview Of Our Terms And Procedures

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## Representative Botanical Reference Materials (rBRMs™)

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## Extract Botanical Reference Materials (XRMs™)

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## Example: RGBRM™ Certificate of Analysis

PRODUCT	<i>Aloe barbadensis</i> (Aloe Vera) Leaves
PART NUMBER	00030822
REFERENCE TYPE	Biomass Reference Material (RG)*
LOT NUMBER	
CDXA NUMBER	
LATIN NAME	<i>Aloe barbadensis</i> P. Mill. [Aloaceae]
COMMON NAME	Aloe Vera
PLANT PART	Leaves
REPORT NUMBER	
DATE OF SAMPLE	07 June 2007
DATE OF REPORT	20 June 2007

\*Note - Reagent Grade (RG) BRM is not vouchered and has not been evaluated by microscopy techniques.

### ANALYTICAL CONDITIONS

TEST	METHOD	RESULT
Macroscopy	Digital Image	Green whole succulent leaves
HP-TLC	CDXA-TLCM-168-00	Conforms (See Below)
Appearance	Visual (Freeze-dried and Milled)	Green fluffy powder

### STORAGE CONDITIONS

**STORAGE** 20-30 °C; Dry storage area; Insect free; Volatile free  
**EXPIRATION DATE** 06/2012 under the above conditions.

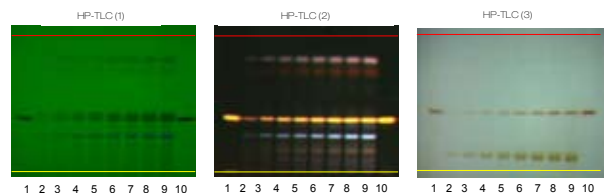
### MACROSCOPY

### HP-TLC CONDITIONS

STATIONARY PHASE	Silica gel 60, F <sub>254</sub> , 10 x 10 cm HP-TLC plates
SAMPLE PREPARATION	0.5 g + 5 ml CH <sub>3</sub> OH sonicate in hot water ~ 10 minutes.
MOBILE PHASE	Ethyl Acetate: Methanol: Water [10/1.4/1]
CHAMBER TEMPERATURE	Ambient
DETECTION	(1) UV 254 nm (2) KOH spray reagent → 365 nm (3) KOH spray reagent → white light

### HP-TLC CHROMATOGRAMS

YELLOW LINE = SAMPLE ORIGIN  
 RED LINE = SOLVENT FRONT @ 60MM

### HP-TLC APPLICATIONS

Lane	ID	Lane	ID
1	Aloin	6	<i>Aloe barbadensis</i> Leaves RGBRM
2	<i>Aloe barbadensis</i> Leaves RGBRM	7	<i>Aloe barbadensis</i> Leaves RGBRM
3	<i>Aloe barbadensis</i> Leaves RGBRM	8	<i>Aloe barbadensis</i> Leaves RGBRM
4	<i>Aloe barbadensis</i> Leaves RGBRM	9	<i>Aloe barbadensis</i> Leaves RGBRM
5	<i>Aloe barbadensis</i> Leaves RGBRM	10	Aloin

Copyright ©2008 ChromaDex™, Inc. All rights reserved.